


InstantBlue®

Protein gel staining in 15 minutes

Get results in just 15 minutes with our ready-to-use InstantBlue® Coomassie protein gel stain.

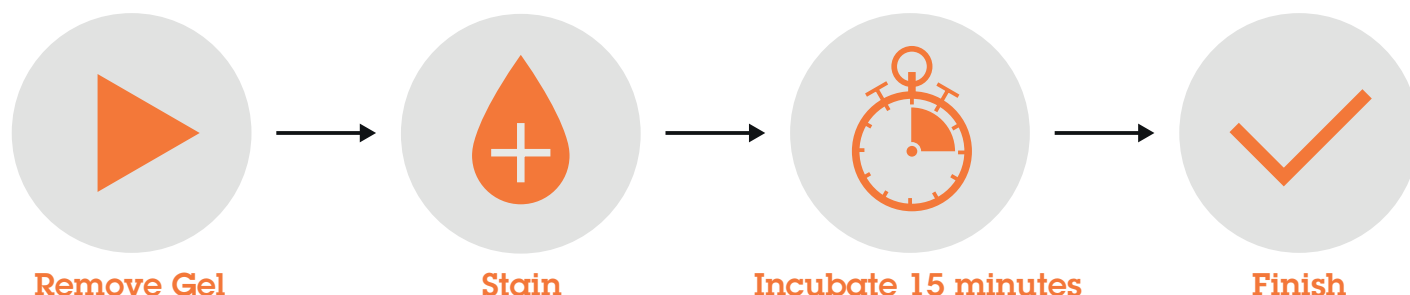

Only proteins are stained, leaving well-defined blue bands on a transparent background with no need to destain and a better signal-to-noise ratio. Gels can remain in the stain for weeks – without the risk of over-staining.

Get back valuable lab time with protein gel staining in 15 minutes.

Benefits of InstantBlue®

- **Ultra-fast staining:** visible bands in 15 minutes or less so you can move your project forward faster
- **Quick and easy to use:** one-step protocol with no washing, fixing, heating, or destaining steps
- **High sensitivity:** detect as little as 5 ng per band of protein (BSA) with clear background
- **Non-toxic formulation:** methanol and acetic acid-free formulation means no gel shrinkage, mass spectrometry compatible, and easy to dispose of

Comparison with other Coomassie stains:

	InstantBlue® (ab119211)	Leading competitor 1	Leading competitor 2	Home-made
No need to wash before staining	✓	3 x 5 minutes	3 x 5 minutes	3 x 5 minutes
No need to fix before staining	✓	✓	✗	✗
No need to heat	✓	✗	✓	✓
No need to destain	✓	✗	✗	✗
No risk of over-staining if left too long	✓	✗	✗	✗
Non-toxic and easily disposed of	✓	✓	✗	✗

Get started with InstantBlue® today.

Contact your sales representative at sales@abcam.com, or scan to learn more:



progress happens together
abcam